

Pyrolysis-GC-HRMS with Soft Ionization


Rachel Sanig

Senior Scientist, *Waters Corporation*

- Introduction to pyrolysis
- Pyrolysis-GC/HRMS with soft ionization
- Applications
 - Comparison to Py-GC/MS with EI
 - Analysis of biobased plastics
 - Statistical analysis
 - Other applications


Combustion

- Hydrocarbon reacts with oxygen to produce carbon dioxide and water

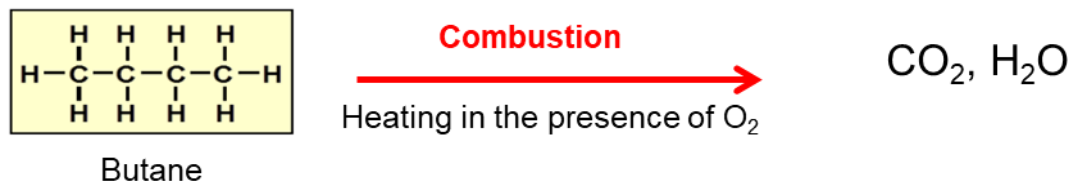

Pyrolysis


- Pyrolysis process is done in the absence of oxygen.

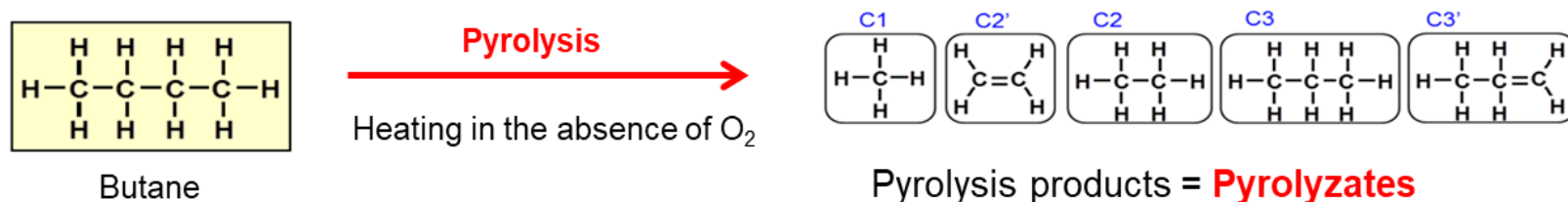

Applied vs Analytical

- There are two key applications for this technique.

Biomass decomposition

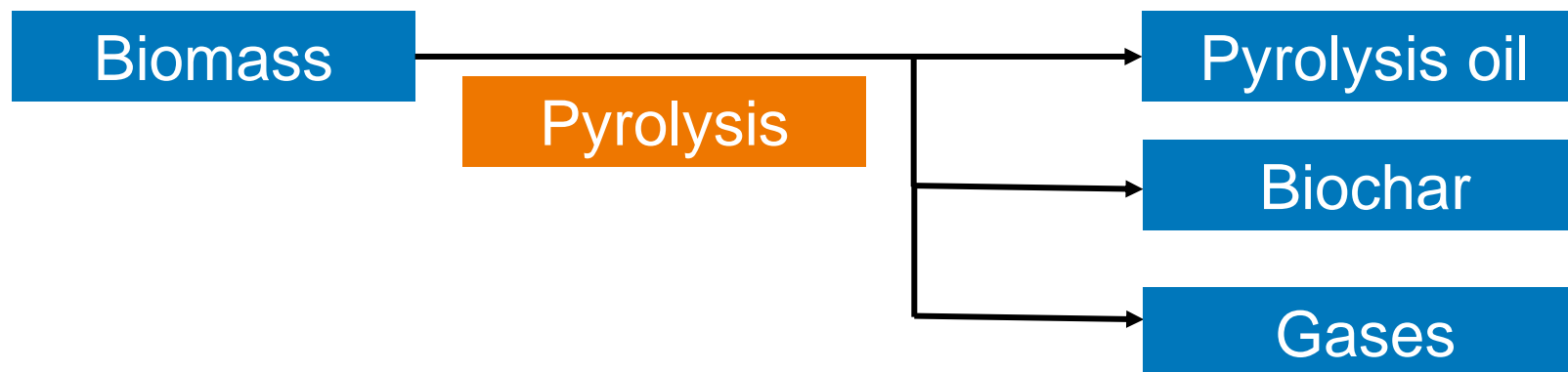


Chemical recycling

- Pyrolysis is **one of the main methods for chemical recycling of plastics**.
 - Breaks downs polymers to monomers with quality recovery of the recycled material
 - Suitable for challenging feedstocks containing mixture of several plastics and impurities

Pyrolysis


- Pyrolysis process is done in the absence of oxygen.

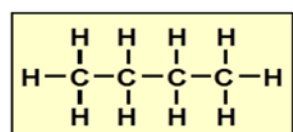

Butane

Pyrolysis

Heating in the absence of O₂


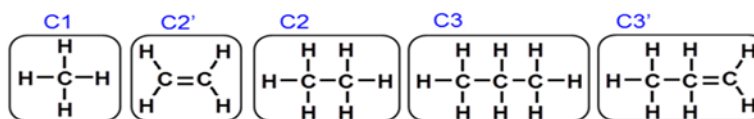


Pyrolysis products = **Pyrolyzates**


GC Separation




Mass Detection

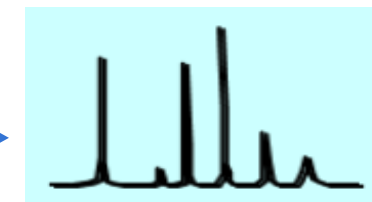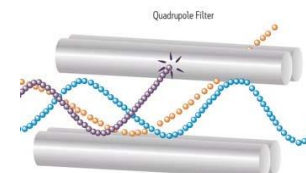


Pyrogram

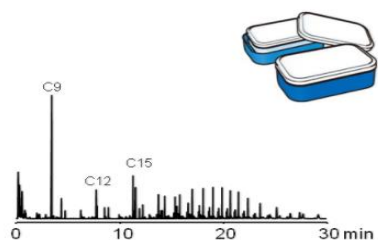

Polypropylene (PP)

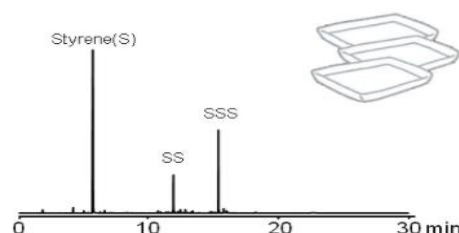

Polystyrene (PS)

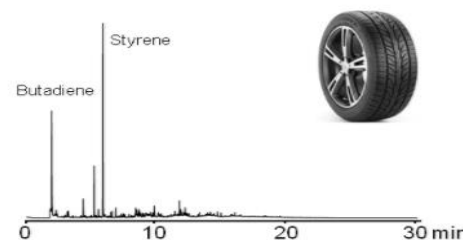

Styrene-butadiene rubber (SBR)

Pyrolysis-GCMS - Challenges and Limitations

■ Database search

- While some monomers are listed in the database, not all pyrolyzates spectrum are registered.

■ Analyzing unknowns

- The pyrolysis process and electron ionization, or electron impact, (EI) do not generate intact molecular ions of pyrolyzates
- Quad mass resolution is also too low to do elucidation and identify true unknowns

■ Sensitivity of high mass range

- High mass pyrolyzates, trimers and higher, are often sensitivity challenged
- EI can make it difficult to confirm the molecular ion of high molecular weight polymers as these often are fragmented in the ionization process

**Py-APGC HRMS
can help with
these analytical
challenges**

Connecting up the instruments




Xevo G2-XS QToF

GC

APGC Source

**CDS Pyroprobe 5000 Series
with autosampler**


**Frontier Auto-Shot Sampler
(AS-1020E)**

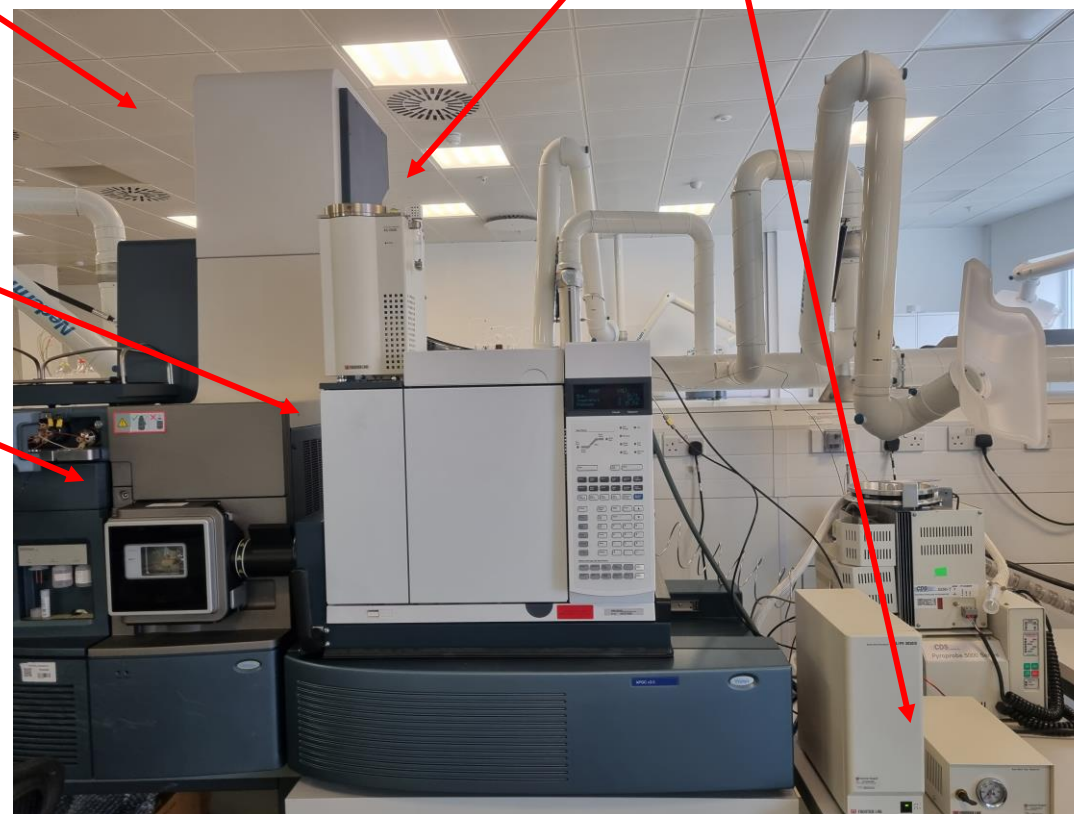

Pyrolysis GC-MS and Soft Ionization

APGC (Atmospheric Pressure Gas Chromatography)

- Soft ionization


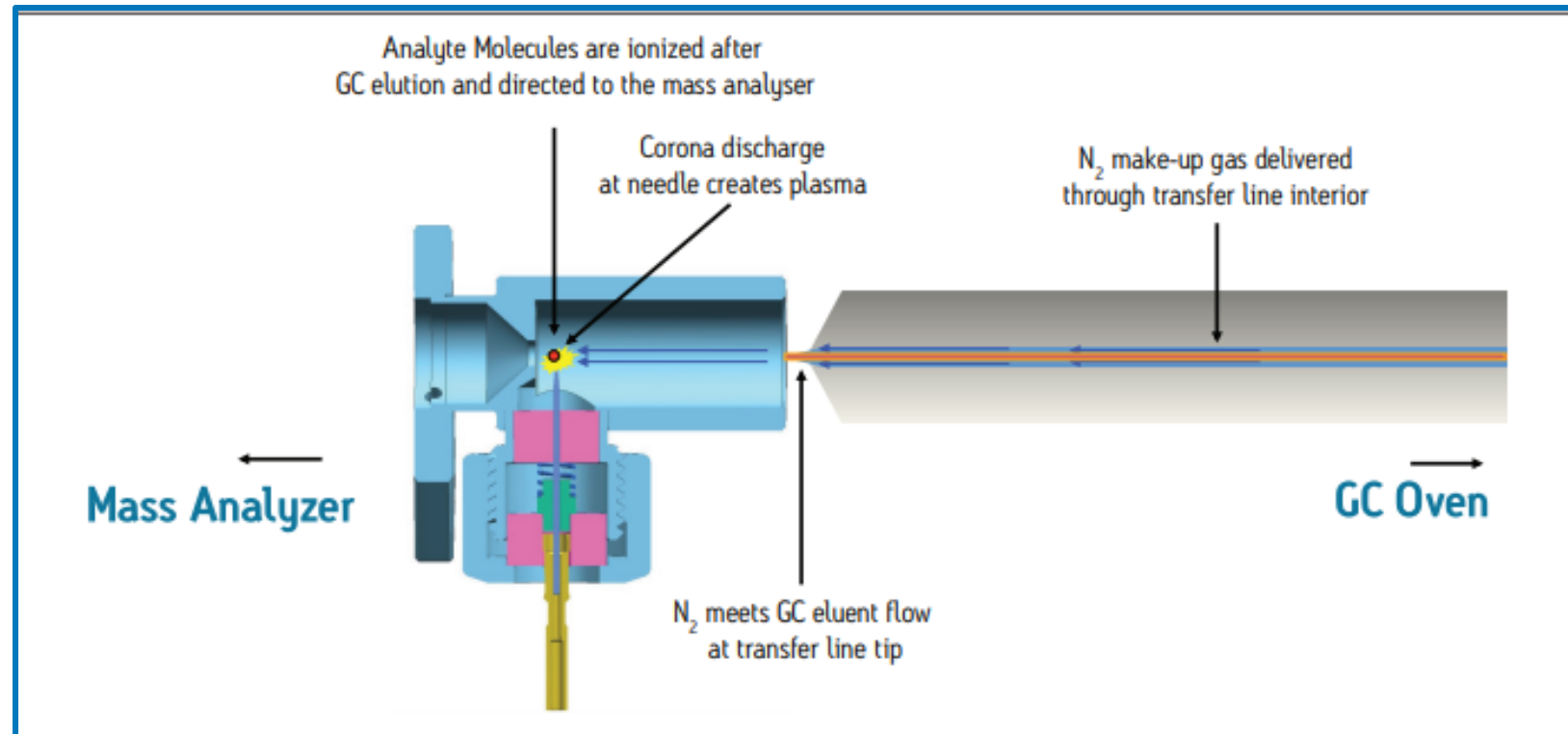


Pyrolysis GC-MS and Soft Ionization

APGC (Atmospheric Pressure Gas Chromatography)

- Soft ionization


“Dry” Source Conditions:
Favored by relatively
non-polar compounds

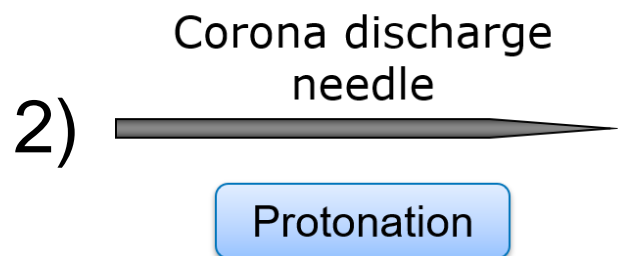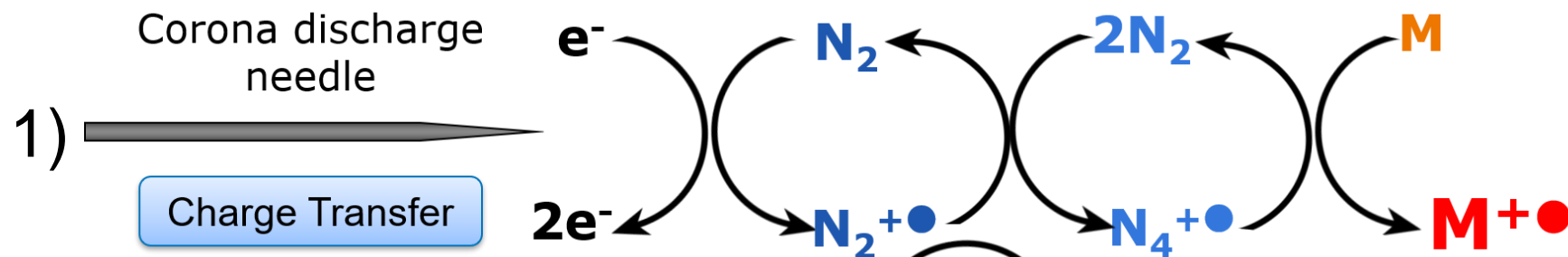

“Modified” Source Conditions:
Favored by relatively
polar compounds

Soft Ionization


- Higher masses can be seen:

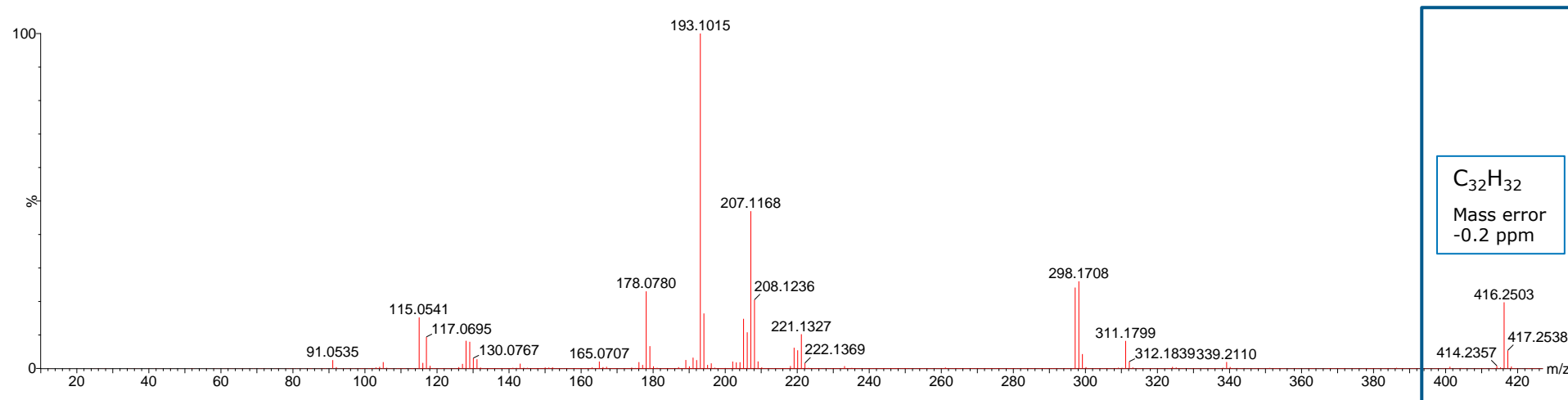

Polystyrene tetramer at m/z 416.2503 corresponding to $[M+C_{32}H_{32}]^+$ cation

Soft Ionization


- Presence of the molecular ion:

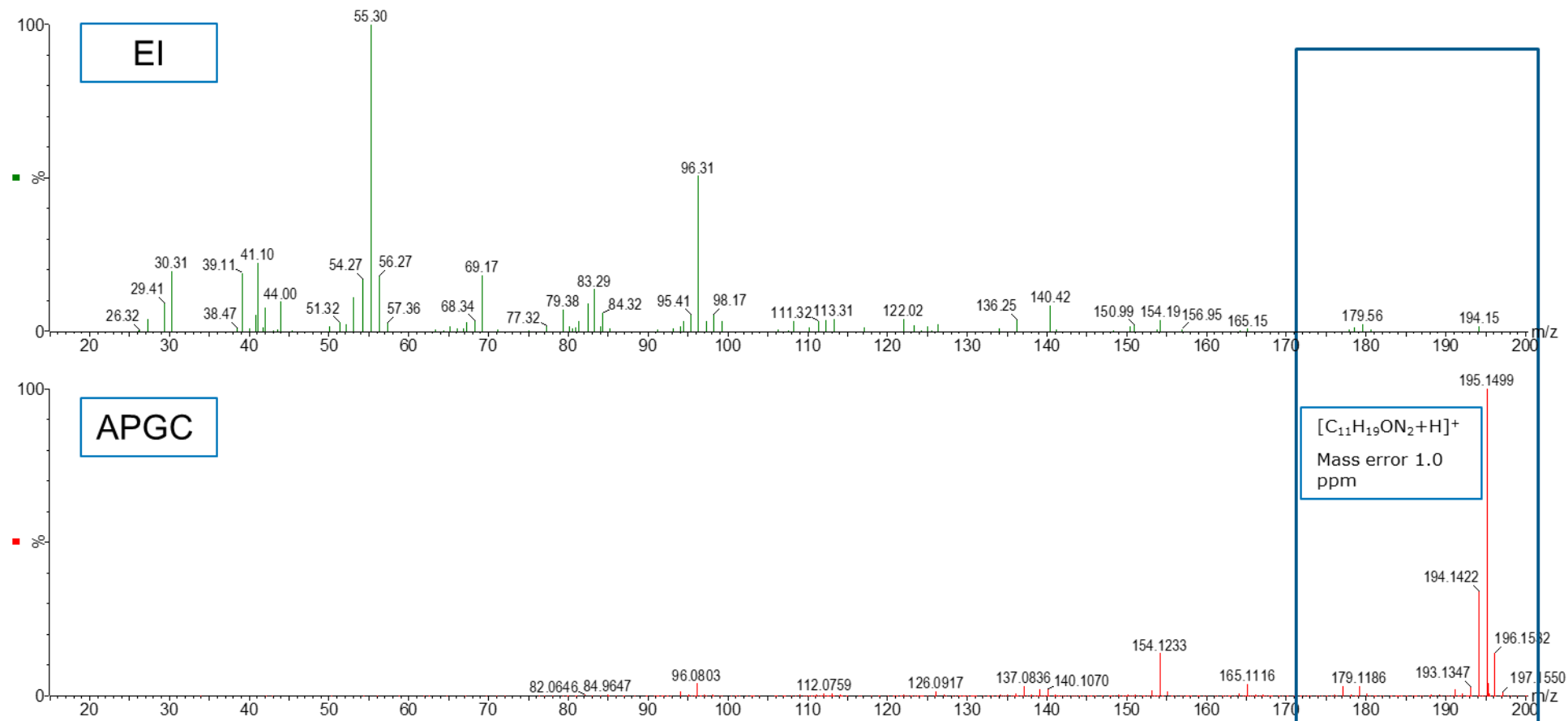

Comparison of spectra from EI and APGC highlighting the presence of the protonated molecular ion in APGC. (polycaproamide (Nylon 6))


MS^E

- High and low collision energies are alternated so that the accurate mass of both precursor and fragments ion can be simultaneously acquired.

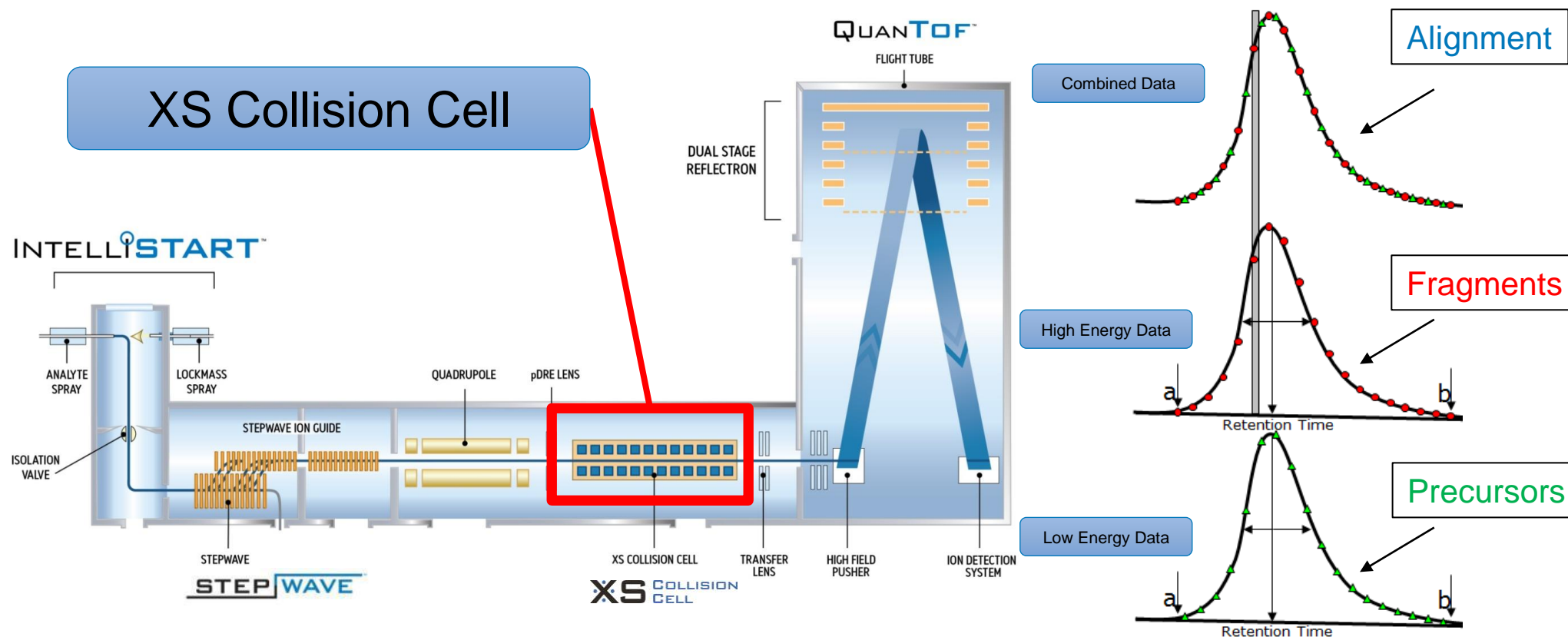


MS^E

- High and low collision energies are alternated so that the accurate mass of both precursor and fragments ion can be simultaneously acquired.


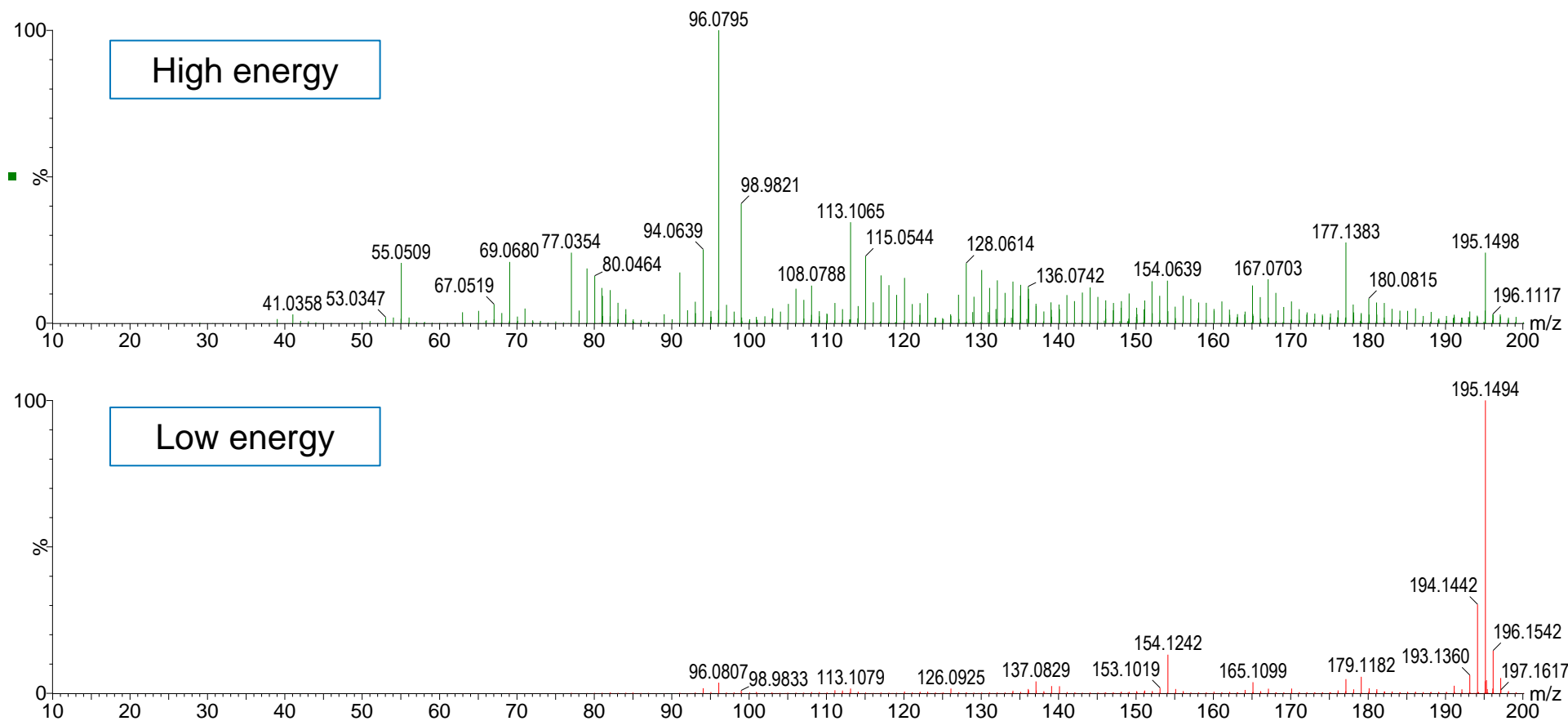


The high and low energy spectra for the protonated ion of m/z 194.2374 using MS^E.

Characterization

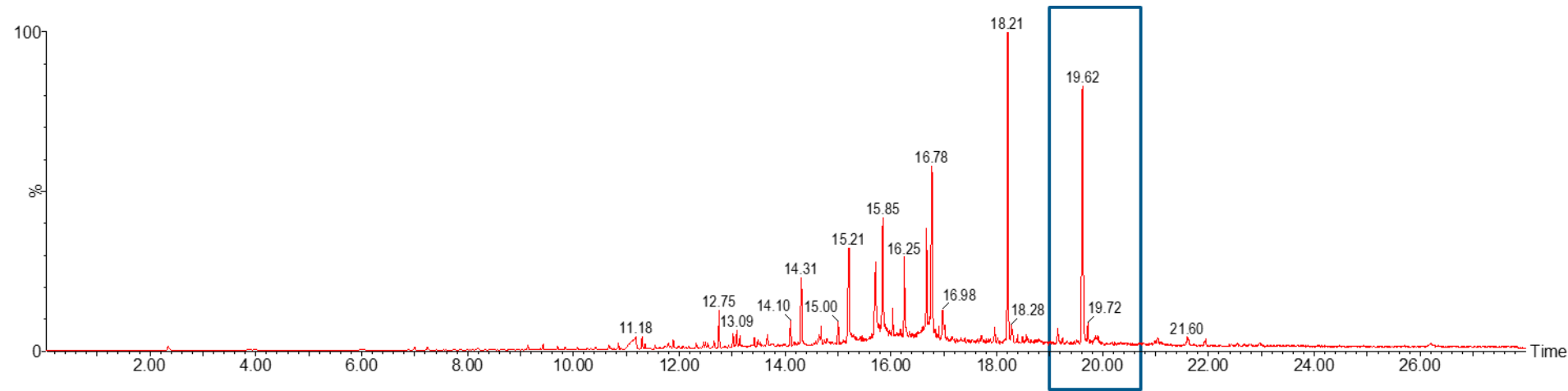




Characterization

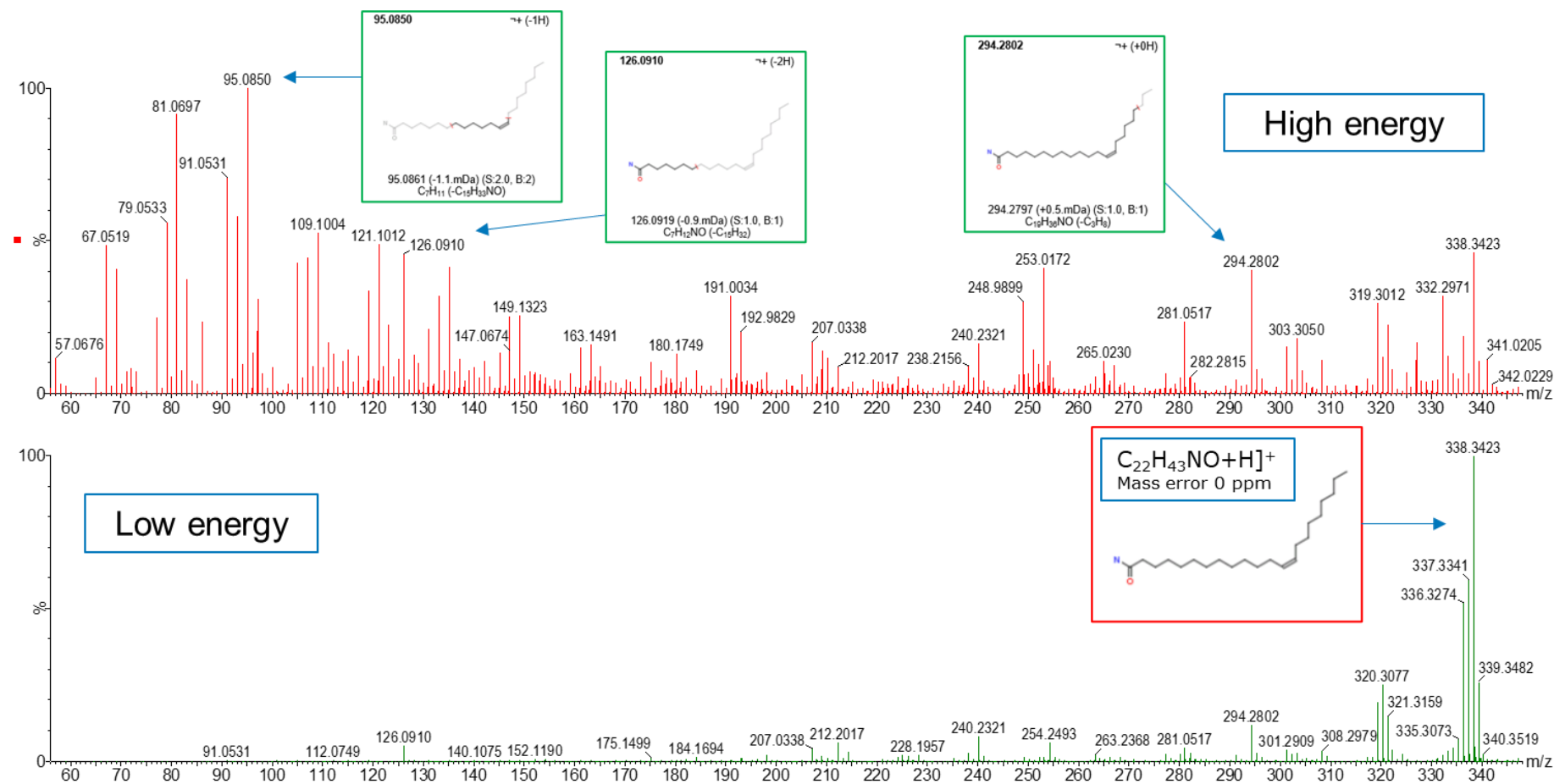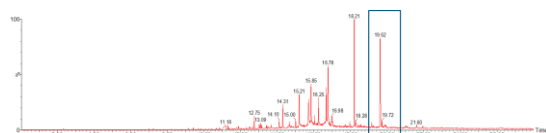

High and low energy spectra for m/z 337.5829 corresponding to erucamide protonated ion using MS^E. Mass fragments assigned by MassFragment.

Characterization of Engineering plastic (PEEK)




- Wet and dry source conditions to understand the polarity of the PEEK pyrolyzates
- Elucidation of polar end groups and components

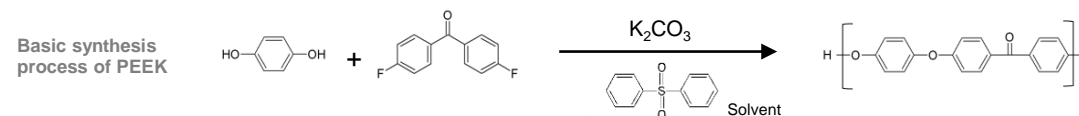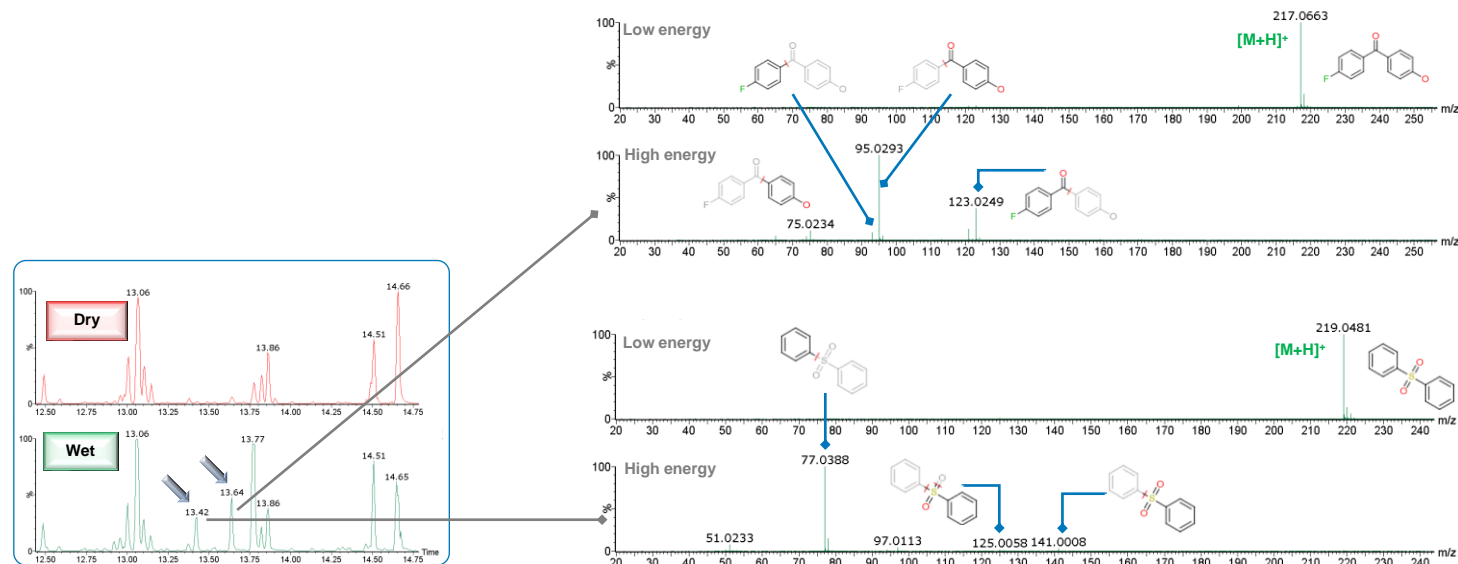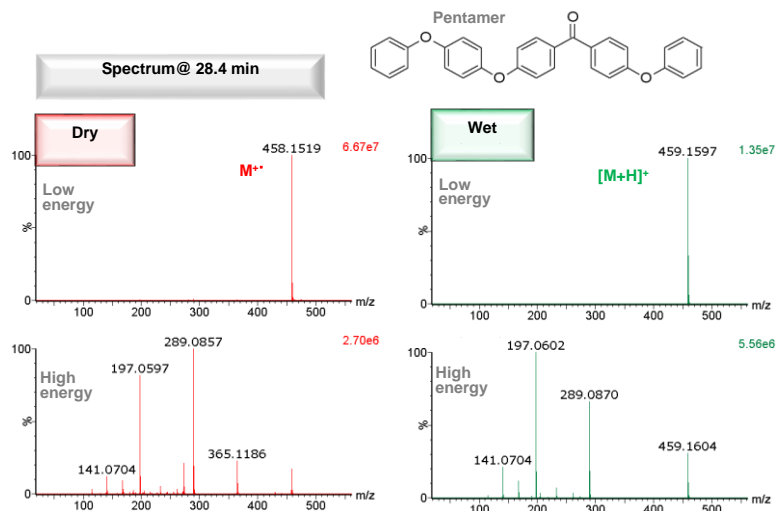

Applications

Comparison of Block and Random Copolymer


- Specific pyrolyzates (markers) were extracted with multivariate analysis

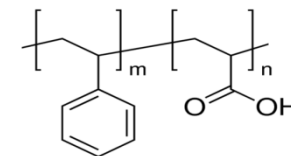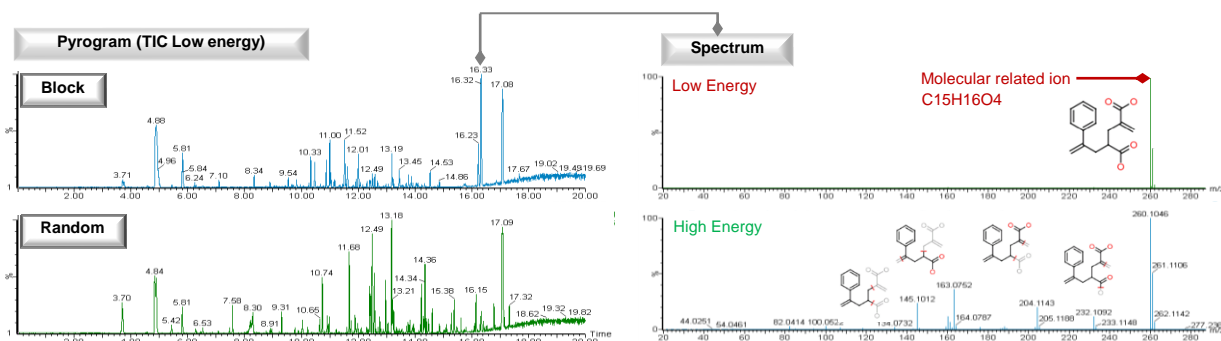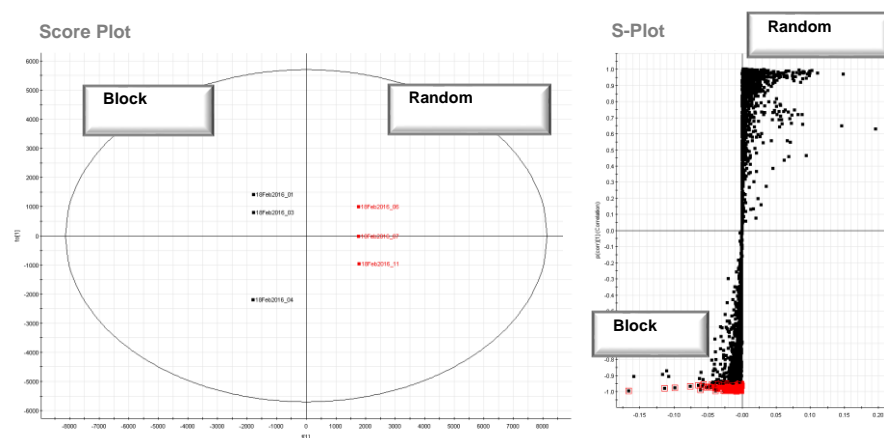


Styrene (Sty) Acrylic Acid (AA)

- Elucidation of pyrolyzates to understand polymer connectivity

Discovered and elucidated markers

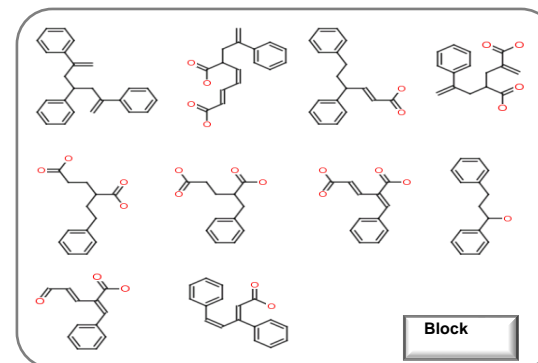

— Sty — Sty — Sty — Sty — Sty — Sty — AA — AA — AA — AA — AA —

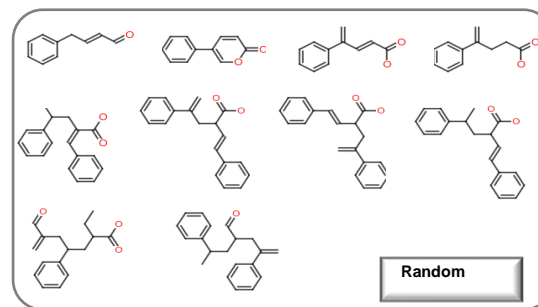

— Sty — AA — Sty — AA — Sty — Sty — AA — Sty — AA — AA — Sty —

Advantages of the Approach

- Analytical pyrolysis-GC-MS for the characterization of complex materials
 - Ongoing project: Recycled materials
- Soft ionization
 - Higher masses
 - Presence of the molecular ion assisting with the identification of unknowns
- MS^E
 - Collection of both the accurate mass of precursor and fragment ions which is key for the structural elucidation of unknowns within a sample. (all of the data, all of the time)
- In house library search/development in addition to existing commercial libraries for putative compound ID
(Poster - TP 138 Combined approach for the characterization of plastics using spectral libraries created from both pyrolysis-GC-MS and pyrolysis-APGC-ToF-MS)